

## Moistening of the Nutrient Pad Sets (NPS)

There are many possible ways of moistening the NPS with the ideal volume of liquid (3.0 – 3,5 ml). We can offer two of these in our product range.

The quick and easy solution are **ampoules** filled with 3.5 ml of sterile, deionised water (article number 6105). Simply twist the top off and pour the content onto the NPS.

The volume of the **self-filling syringe** (article number 6100) can be infinitely varied up to 5.0 ml. The pre-set volume is exactly dosed sterile onto the NPS through a sterile syringe filter unit with a pore size of 0.2 µm. The syringe is then re-filled automatically with deionised or demineralised water from a reservoir.

Please contact us in case of any questions. We will be happy to assist you.

**Dr. Möller& Schmelz GmbH**  
**Gesellschaft für angewandte Mikrobiologie**

Robert-Bosch-Breite 15  
D-37079 Göttingen

Germany

☎ +49 (0)551 66708

☎ +49 (0)551 68895

✉ info@moeller-schmelz.de

www.moeller-schmelz.de